

Priming Procedure

Startup Guide

Preparing Your Pump for First Time Use




Self-Flush

- Connect solvent inlet and outlet tubing as shown.
- Attach syringe to 1" (clear urethane) tubing
- Draw syringe back to prime.
- Remove syringe and place tubing in solvent. *Replace solvent weekly*

Note: If self-flush option is not used, refer to manual for proper pump modification.

Pump

- Connect pump inlet tubing.
Make sure ferrule is in the correct position
- Turn Power on
- Attach syringe to Prime / Purge.
- Open Prime / Purge valve by turning knob counterclockwise two turns.
- Draw syringe back to prime
Draw approximately 15 mL of fluid.
- Press PRIME button, continue to draw on syringe until no bubbles are seen,
- Close Prime / Purge valve
- Press PRIME button
- Remove syringe
Replace solvent weekly

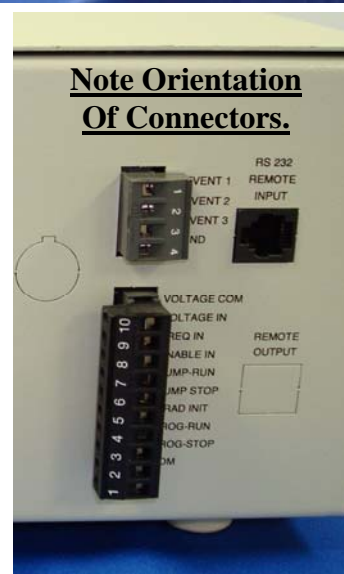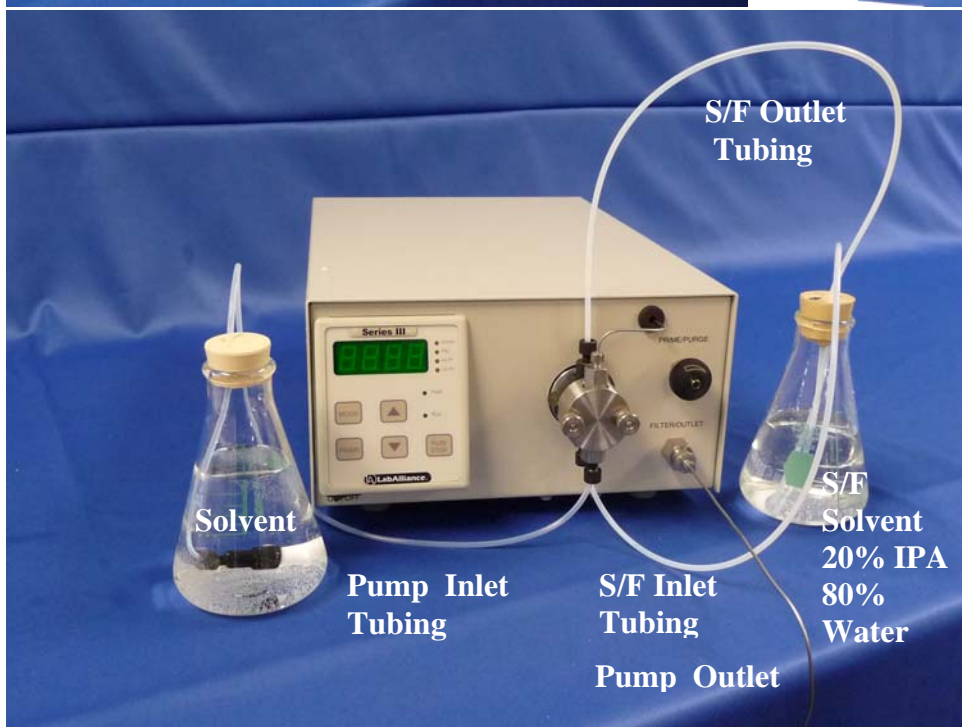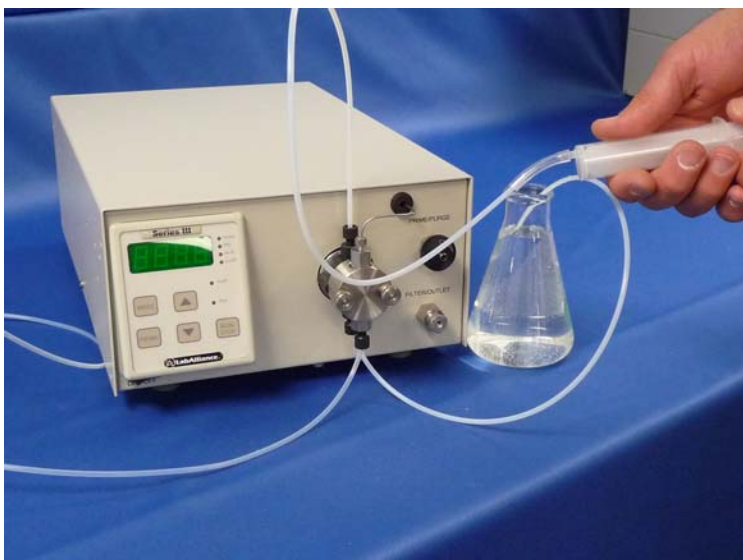

Note Orientation Of Connectors.